

Remove the PS5000 Slide Jam SOP

You can find the step-by-step video on our website

<http://www.scanace.com>

1. Lift GUIDE PLATE to find the jammed slide.




2. Use a small tweezer type tool to grab the jammed slide and push it to the slide scan window. Maneuvering it in small increments back into the scanner



3. Push the slide until it is at the scan window proper position.



4. Push the Green button on top to open the Single Slide Release to raise jammed slide and take it out.

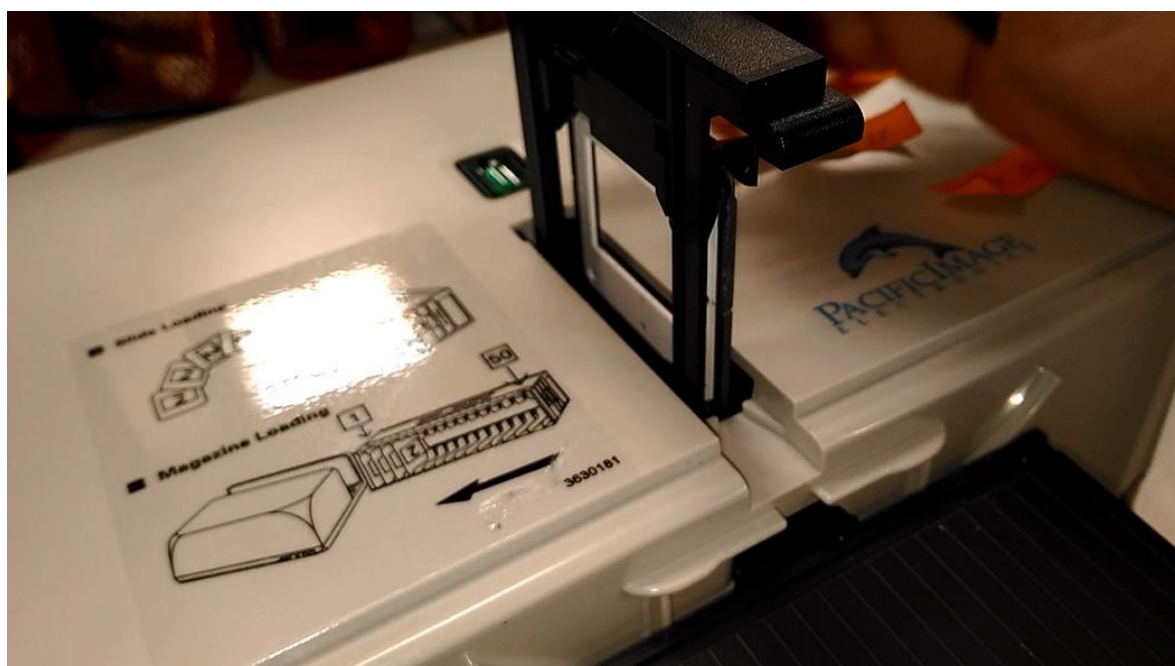

5. Pull the slide transport arm out, and take out the Slide Tray to resolve the slide jam issue.

